Installation

S19-899

On-Site® Gravity-Fed Eye Wash Refill Kit

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Installation

Fill and mount tank

1. Pour clean, potable water into the tank; continue pouring until the tank is 1/3 to 1/2 full.
2. Pour the entire contents of the water preservative bottle into the eye wash tank.
3. Add clean, potable water to the tank until the water level reaches the top of the tank.
4. Insert one liner into the cap with hole and attach the cap to the top of the eye wash tank (make sure the cap is tightly closed) (see Figure 1).

- S19-881 ONLY: Insert the tube into the pinch to prevent leaking when installing the tank.

NOTE FOR TANK INSTALLATION: To ensure proper installation, the tank should be placed on the pedestal from the front of the unit.

5. Invert the tank and place it on top of the pedestal as shown in Figure 2a. The pierce on the pedestal will puncture a hole in the tank cap liner and allow the solution to flow.
 - S19-881 ONLY: Make sure that the locking tabs are in front of the pedestal (see Figure 1). Raise the arm to snap fit in an upright position (this will automatically release the pinch).
 - S19-921 ONLY: Raise the arm to snap fit in an upright position, then secure the tank to the pedestal using the side tank clamps. Make sure the latches are pulled down tight to the pedestal (see Figure 2b).

NOTE: After the tank is mounted to the pedestal, minor spillage could occur. Allow the fluid to stabilize completely before inspecting the unit for leaks.

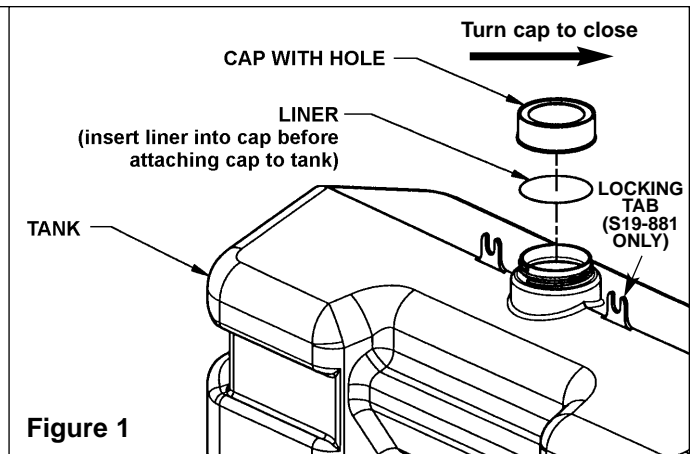
Figure 1

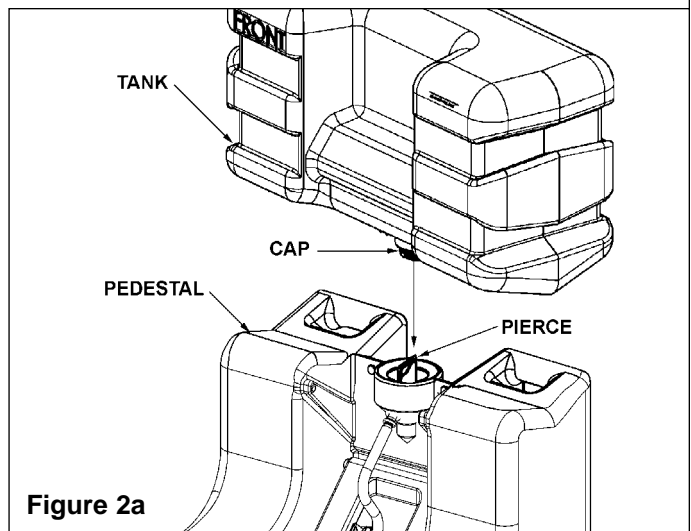
Figure 2a

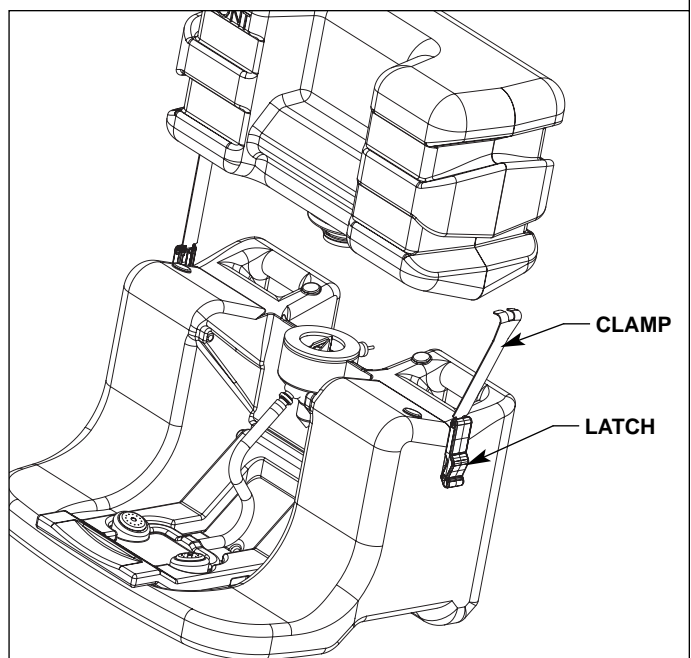
Figure 2b

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